



Sartobran® P 0.1 µm

Sterilizing Grade Filter Cartridges



Description

Sartobran® P 0.1 µm rated, high-flow filter elements are designed to give enhanced sterility assurance for applications with microorganisms present that can pass through 0.2 µm rated sterilizing grade filters. The Sartobran P's cellulose acetate membrane offers ultra low binding properties for proteins and preservatives, making Sartobran P filters the ideal choice for filtration of high value bio-pharmaceutical products.

Applications

All applications which require sterilizing grade filtration with retention finer than conventional 0.2 µm sterilizing grade filters for removal of unusually small microorganisms. This typically includes:

- Bio-processed pharmaceuticals
- Long term filling operations
- Filtration in pharmaceutical water systems

Flow Rates

Higher flow rates than other 0.1 µm rated filters provide short filtration time and gentle product treatment, even if replacement of conventional 0.2 µm rated filters is necessary.

Total Throughput

Due to the "built-in prefiltration" by a 0.45 µm membrane, Sartobran P 0.1 µm filters provide higher total throughputs than any other 0.1 µm rated filter for economical process design.

Highest Product Yield

The ultra low adsorption characteristic of the Sartobran P's cellulose acetate membrane provides the highest product yield – especially important for high value proteins.

Flexibility

Sartobran P 0.1 µm filters are available in traditional cartridge formats and disposable capsules from 500 cm² to 1.8 m² for simple linear scale up and process flexibility.

Quality Control

Each individual element is integrity tested by diffusion and bubble point test prior to release, assuring absolute reliability.

Documentation

Sartobran P cartridges are designed, developed and manufactured in accordance with an ISO 9001 certified Quality Management System. A Validation Guide and Extractables Guide are available for compliance with regulatory requirements.

Specifications

Materials

Prefilter Membrane:	Cellulose Acetate
Endfilter Membrane:	Cellulose Acetate
Support Fleece:	Polypropylene
Core:	Polypropylene
End Caps:	Polypropylene
O-Rings:	Silicone (optional EPDM or Viton)

Pore Size

0.45 µm + 0.1 µm

Available Sizes | Filtration Area

Cartridges

Size 1	10"	0.6 m ²	6.5 ft ²
Size 2	20"	1.2 m ²	12.9 ft ²
Size 3	30"	1.8 m ²	19.4 ft ²

Mini Cartridges

Size 7	0.05 m ²	0.54 ft ²
Size 8	0.1 m ²	1.1 ft ²
Size 9	0.2 m ²	2.2 ft ²

Available Adapters Cartridges

21, 25, 27, 28

Available Adapter Mini Cartridges

15

Operating Parameters

Max. Allowable Differential Pressure:	5 bar 75 psi at 20 °C
	2 bar 29 psi at 80 °C
Max. Allowable Back Pressure:	2 bar 29 psi at 20 °C

Specifications

Extractables

Sartobran® P 0.1 µm rated filter cartridges meet or exceed the requirements for WFI quality standards set by the current USP.

Regulatory Compliance

100% Individually integrity tested

Integrity test correlated to HIMA/ASTM F 838-05 Bacteria Challenge Test.

Non-pyrogenic according to USP Bacterial Endotoxins

Passes USP Plastics Class VI Test

Non-fiber releasing according to 21 CFR

Sterilization

In-Line Steam Sterilization

134 °C, 20 min. at max differential pressure of 0.5 bar | 7 psi

Autoclaving

134 °C, 2 bar | 29 psi, 30 min

Sterilization Cycles

In-Line Sterilization: Min. 25
(only cartridges)
Autoclaving: Min. 25

Technical References

Validation Guide:

SPK5726-e


Extractables Guide:

SPK5720-e

Ordering Information


Order Code	Pore Size [µm]	Pack Size [Pieces]	Test Pressure [bar psi]	Max. Diffusion [ml/min]	Min. Bubble Point [bar psi]
Standard Cartridges					
523**58H1----P	0.1	1	3.0 44	15	3.8 55
523**58H2----P	0.1	1	3.0 44	30	3.8 55
523**58H3----P	0.1	1	3.0 44	45	3.8 55
Mini Cartridges					
5231558H7----B	0.1	5	3.0 44	3	3.8 55
5231558H8----B	0.1	5	3.0 44	4	3.8 55
5231558H9----B	0.1	5	3.0 44	5	3.8 55

Water Flow Rates for 10", 20" and 30" Cartridges

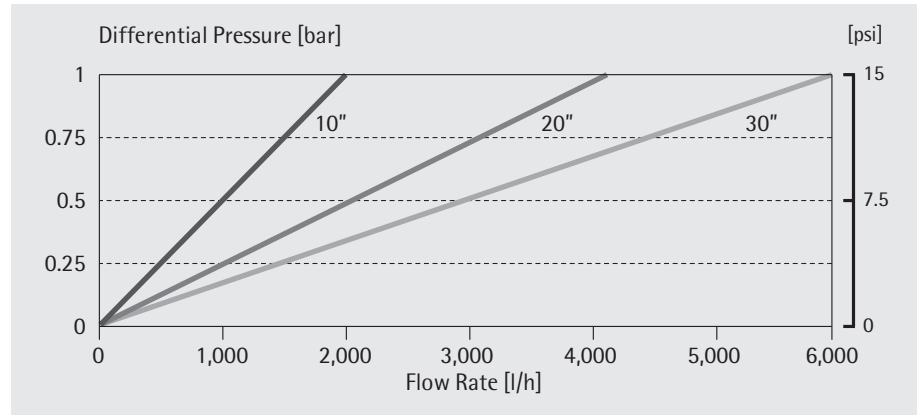

Standardized at 20 °C

Total Throughput Comparison

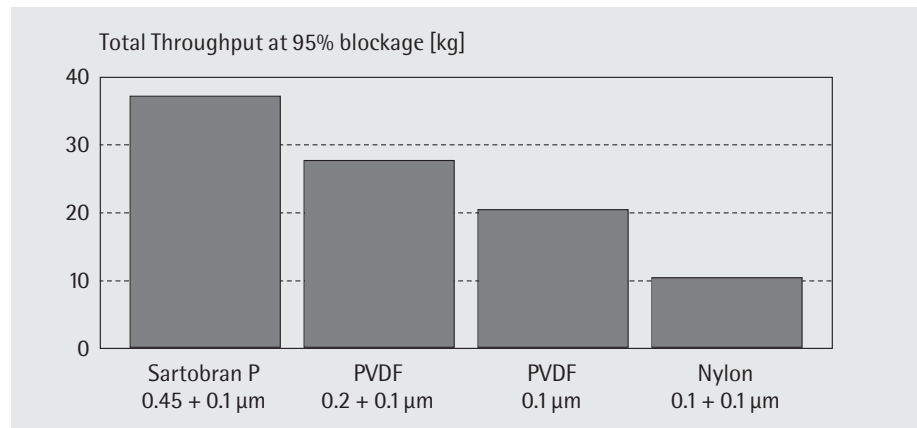

10" Cartridge format

Sartorius Stedim Biotech GmbH
August-Spindler-Strasse 11
37079 Goettingen, Germany

Phone +49.551.308.0
Fax +49.551.308.3289
www.sartorius-stedim.com

USA Toll-Free +1.800.368.7178
UK +44.1372.737159
France +33.442.845600
Italy +39.055.63.40.41
Spain +34.91.3586102
Japan +81.3.3740.5407

Specifications subject to change without notice. Printed and copyrighted by Sartorius Stedim Biotech GmbH W · G
Publication No.: SPK2041-e09044
Order No.: 85030-513-73
Ver. 04 | 2009